

Instructions for sending molecular materials for deposit

Patent deposit

1. Deposit form

Submit a signed deposit form together with the molecular material.

2. Molecular material

2.1. Molecular materials are contained in cryotubes ready for preservation by freezing.

2.2. Molecular material name and date of preservation must be printed on self-adhesive stickers using laser printer ink. Each label should be further secured with a layer of sealing film or clear tape.

2.3. During transportation, molecular material must be maintained at -70°C to -80°C.

2.4. Send 11 tubes per sample.

2.5. Molecular materials classified in Risk Group 2 must be packed and shipped according to national biosafety and biosecurity regulations or IATA Dangerous Goods Regulations.